Mikroskop · Microscope · Microscopio

## **Biorit TP**

Art. No. 5101100

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**EN** Instruction Manual

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<b>Deutsch .....</b>	<b>3</b>
<b>English.....</b>	<b>14</b>

# 1 Impressum

Bresser GmbH

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46414 Rhede

Deutschland

<http://www.bresser.de>

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Deutschland

<http://www.bresser.de>

For information regarding liability claims or service requests, please refer to chapters "Warranty" and "Service" in this documentation. We ask for your understanding, that any requests or submissions that have been addressed to the manufacturer's address can not be processed.

Errors reserved - technical specifications subject to change.

Copyright: 2016Bresser GmbH

All rights reserved.

It is not allowed to reproduce this documentation or parts of it in any kind (e.g. photocopying, printing, etc) without written a permission of the manufacturer. It is also prohibited to use or transmit this documentation with electronic systems (e.g. image file, website, etc).

The descriptions and brand names of the corresponding firms used in this documentation are generally protected under trademark, brand and/or patent law for Germany, for the European Union and/or other countries.

---

<b>1</b>	<b>Imprint</b>	<b>14</b>
<b>2</b>	<b>Features</b>	<b>16</b>
<b>3</b>	<b>About this Instruction Manual</b>	<b>16</b>
<b>4</b>	<b>Intended Use</b>	<b>16</b>
<b>5</b>	<b>General safety instructions</b>	<b>16</b>
<b>6</b>	<b>Product images</b>	<b>18</b>
<b>7</b>	<b>Scope of delivery</b>	<b>18</b>
<b>8</b>	<b>Parts overview</b>	<b>18</b>
<b>9</b>	<b>Site location</b>	<b>19</b>
<b>10</b>	<b>Installing the USB power adapter</b>	<b>19</b>
<b>11</b>	<b>Setting up power supply</b>	<b>20</b>
<b>12</b>	<b>Optional power supply</b>	<b>20</b>
<b>13</b>	<b>Turning on/off the device</b>	<b>20</b>
<b>14</b>	<b>Adjusting the lighting</b>	<b>21</b>
<b>15</b>	<b>Inserting slides</b>	<b>21</b>
<b>16</b>	<b>Observation</b>	<b>22</b>
<b>17</b>	<b>Using the condenser</b>	<b>22</b>
<b>18</b>	<b>Using the iris diaphragm</b>	<b>23</b>
<b>19</b>	<b>Cleaning and maintenance</b>	<b>23</b>
<b>20</b>	<b>Disposal</b>	<b>23</b>
<b>21</b>	<b>Technical data</b>	<b>24</b>
<b>22</b>	<b>Warranty</b>	<b>24</b>
<b>23</b>	<b>EC Declaration of Conformity</b>	<b>24</b>

## 2 Features

- Biological microscope for transmitted light microscopy
- Optional off-grid energy supply with integrated battery
- Cross table with Nonius scale and with lateral and longitudinal adjustment
- LED illumination with stepless brightness adjustment
- Objectives and eyepieces exchangeable
- Condensor with filter holder
- Universal USB power adapter with two adapter attachments

## 3 About this Instruction Manual

### NOTICE

**These operating instructions are to be considered a component of the device.**

Please read the safety instructions and the operating instructions carefully before use.

Keep these instructions for renewed use at a later date. When the device is sold or given to someone else, the instruction manual must be provided to the new owner/user of the product.

## 4 Intended Use

- This device is not intended for use by individuals (including children) with limited physical, sensory or mental capabilities or those lacking in experience and/or knowledge, unless they are supervised by an individual responsible for their safety or have received instructions from them regarding the use of the device.
- This product is intended only for private use.
- It was developed for the magnified display of things in nature.
- The device is intended only for indoor use.

## 5 General safety instructions

### DANGER

**Danger of an electric shock!**

This device contains electronic components that operate via a power source (power supply and/or batteries). In case of any improper use of this device, there is a risk of an electric shock. An electric shock can cause severe injury or even death. Therefore please read the safety instructions below to avoid an explosion.

- Disconnect the device from the power supply by pulling the power plug when it is not used or in case of longer interruption of operation and before starting any work on maintenance and cleaning.
- Position your device so that it can be disconnected from the power supply at any time. The wall socket you use should be located near the device and easily accessible, since the plug on the power cable serves as a disconnecting device for the power supply.
- Always pull on the plug to separate the device from the power supply. Never pull on the cable.
- Before operating, check the device, cables and connections for damage.
- Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.

- Only use the device in complete dry environment and do not touch it with wet or moist parts of your body.

## **DANGER**

### **Danger of choking!**

In case of any improper use of this device, there is a risk of choking, especially for children. Therefore please read the safety instructions below.

- Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard!

## **DANGER**

### **Danger of explosion!**

In case of any improper use of this device, there is a risk of an explosion. Therefore please read the safety instructions below to avoid an explosion.

- Do not expose the device to high temperatures. Use only the supplied power adapter. Do not short-circuit the device or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.

## **CAUTION**

### **Danger of injury!**

This device contains components and/or accessories that can cause minor to severe injuries in case of any improper use. Therefore please read the safety instructions below to avoid any bodily injury.

- Tools with sharp edges and points are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.
- Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after using the chemicals. In the event that the chemicals come into contact with your eyes or mouth, rinse thoroughly with water. If you are in pain after exposure, contact a doctor immediately and take the substances with which you came into contact with you.

## **CAUTION**

### **Fire hazard!**

In case of any improper use of this device, there is a risk of fire. Therefore please read the safety instructions below to avoid the initiation of burning.

- Never cover the ventilation slots or cooling fins of the device while using it or as long as it has not sufficiently cooled down!

## **NOTICE**

### **Risk of property damage!**

In case of any improper use of this device and/or its accessories, there is a risk of property damage. Therefore only use the device according to the safety instructions below.

- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Do not expose this device to higher temperatures and protect it from water and high humidity.

- Protect the device from severe shocks!
- For this device only use accessories and spare parts that comply with the technical information.

## 6 Product images

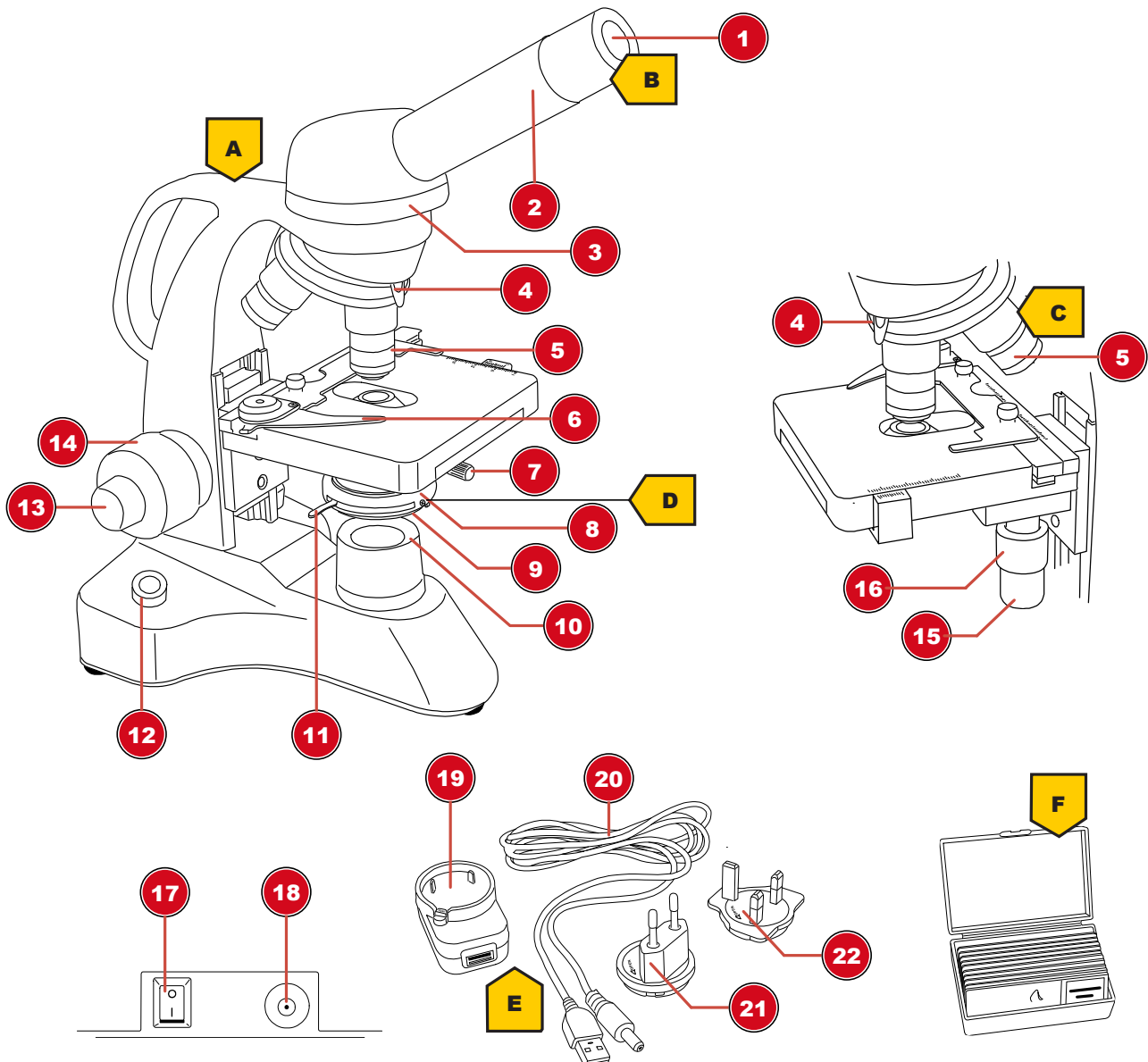
Illustration 1: Product images

## 7 Scope of delivery

Microscope (A); 10x WF Eyepiece (B); 3 Objectives (pre-assembled) (C); Condenser (pre-assembled) (D); Power adapter with plug and cable (E); Box with preparations, blank slides and cover glasses (F)

## 8 Parts overview

1	Eyepiece
2	Eyepiece holder
3	Microscope head
4	Objective nosepiece

5	Objective (4x, 10x or 40x)
6	Specimen clamp
7	Condenser height adjustment
8	Condenser
9	Filter holder
10	Illumination
11	Adjusting lever for iris diaphragm
12	Rotary knob for brightness adjustment
13	Fine focusing
14	Coarse focusing
15	Cross table lateral adjustment (X-axis)
16	Cross table longitudinal adjustment (Y-axis)
17	On/Off switch
18	DC power port
19	Universal USB power adapter
20	USB power cable with coaxial/barrel connector
21	Power adapter attachment (EU)
22	Power adapter attachment (UK)

## 9 Site location

Choose a suitable location before assembly and commissioning.

Place the device on a stable and flat surface, free from vibrations.

## 10 Installing the USB power adapter

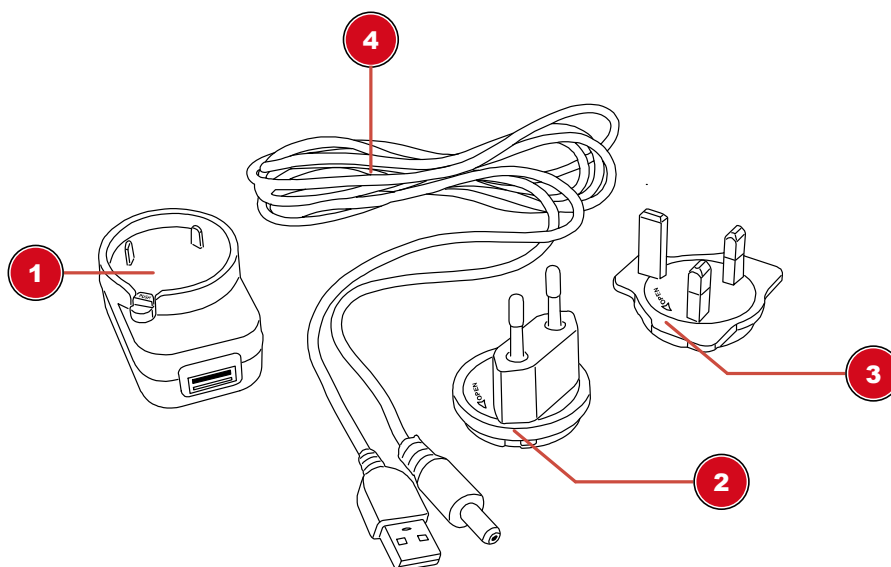
Illustration 2: USB power adapter with exchangeable attachment

1	USB power adapter	2	Adapter attachment (EU)
3	Adapter attachment (UK)	4	USB power cable with coaxial/barrel connector USB power cable with coaxial/barrel connector

1. Place suitable adapter attachment with its notches on the pins of the USB power adapter.
2. Turn adapter attachment in LOCK direction until it is locked.

3. Put the USB plug of the power cable into the USB port of the power adapter.
4. To remove the adapter attachment, push the PUSH button and turn the attachment in OPEN direction.

## NOTICE

### Using the USB power cable with a computer

Alternatively the device can be powered by a computer via the included USB power cable: Put the USB plug into a free USB port of the computer.

## 11 Setting up power supply

Illustration 3: DC connection socket for coaxial/barrel connector

1 On/Off switch

2 DC connection socket

1. Put the DC connector of the power cable into the DC connection socket of the device.
2. Put the mains plug of the power cable into the power outlet.

## 12 Optional power supply

## NOTICE

### Avoid deep discharging of the rechargeable battery!

If the light will be weaker without any reason, it indicates that the battery power is low. Then the the device must be connected to a power supply immediately to avoid deep discharging of the rechargeable battery. Deep discharging can cause irreparable damage to the battery and the electronics.

This device can be operationed alternatively via the integrated battery without need of a power supply. When disconnecting the device from the power supply, it switches to battery powered operation automatically.

## 13 Turning on/off the device

1. Move On/Off switch to position 'I' to turn on the device.
2. Move On/Off switch to 'O' position to turn off the device.

## 14 Adjusting the lighting

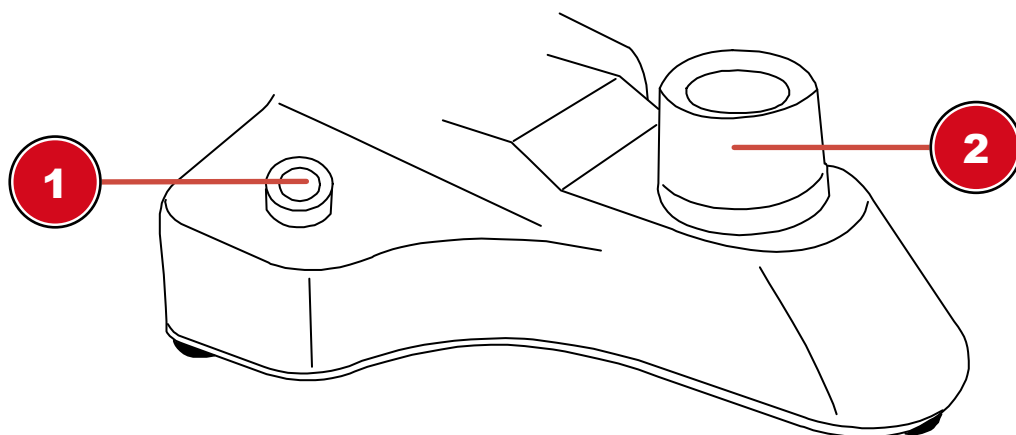
Illustration 4: Transmitted light unit with dimmer

1 Dimmer control

2 Lighting unit

This device provides an illumination with a permanently installed LED light source.

### NOTICE

**Do not operation the device on highest brightness over a longer time!**

This will reduce the lifetime of the LED.

To control the brightness proceed as follows:

1. Turn on the device.
2. Turn/Move the dimmer control to adjust the brightness of the light source.

### NOTICE

**Illuminating the object optimally**

Adjust the brightness of the light source so that the illumination works at its optimum. Depending on light influences from outside and the object being observed, the illumination must be adjusted more often.

## 15 Inserting slides

1. Move the object plate downwards completely by turning the focus wheel.
2. Press the object clamping outwards.
3. Insert the slide straight into the fixture's angle.
4. Let the object clamping fall back carefully until the slide is fixed in the fixture.

## 16 Observation

### NOTICE

#### Using eyepieces and objectives

When using a specific eyepiece or objective, you will get the individual magnification of an object being observed. The magnification is imprinted on the eyepiece and objective with a large number (e.g. 10x for 10-fold magnification, 4 for 4-fold magnification).

1. Move the object plate downwards completely by turning the focus wheel.
2. Insert a slide and place it in the middle under the objective.
3. Set an objective with lowest magnification below the object plate.
4. Insert an eyepiece with lowest magnification in the eyepiece holder.
5. Turn on the device and the illumination.
6. Look through the eyepiece and move the object plate upwards by turning the focus wheel until a sharp image is visible.

### NOTICE

#### Avoid damages to the specimen!

Only turn the focus wheel until it stops to avoid damages to the specimen throughout collision with the objective.

1. If necessary, adjust the image sharpness by turning the fine focus adjustment.

## 17 Using the condenser

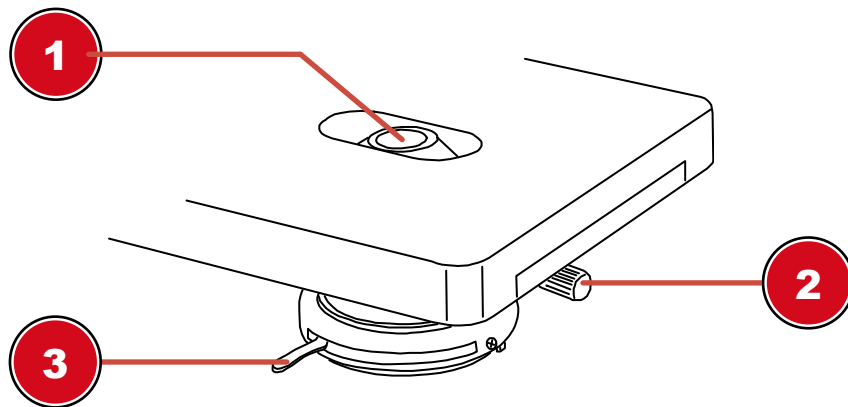
Illustration 5: Condenser

1 Condenser lens

2 Condenser height adjustment

3 Adjusting lever for iris diaphragm

The condenser is used to adjust the light bundling to the observed object in an optimal way. Light bundling can be varied by moving the condenser height adjustment.

## 18 Using the iris diaphragm

### NOTICE

**The iris diaphragm is used to increase the depth of focus.**

With an open aperture more light is getting through but the depth of focus is less. With a partially closed aperture the depth of focus is higher but it is necessary to increase the brightness of the illumination then.

It is recommended not to close the iris diaphragm too much when working with higher magnifications. Otherwise the image resolution will decrease.

Move the lever for the iris diaphragm to right or left to change the aperture setting.

## 19 Cleaning and maintenance

- Protect the device against dust and moisture.
- Avoid putting fingerprints and other contaminations on any optical surfaces.
- Make sure the residual moisture is fully eliminated before use.
- Disconnect the device from the power supply by pulling the power plug when it is not used or in case of longer interruption of operation and before starting any work on maintenance and cleaning.
- Only use a dry cloth to clean the exterior of the device. To avoid damaging the electronics, do not use any cleaning fluid.
- Do not use organic solvents (e.g. alcohol, ether, acetone, xylol or other dilutions) to clean lacquered parts or plastic parts!
- Disassemble optical elements (e.g. objective, eyepiece, etc) before cleaning.
- Blow away loose dust from the lens surfaces first.
- Use high quality lens cleaning towels or soft cloth and moisten it with a little bit of pure alcohol (available in a drug store). Wipe the lens surface with it.
- Pull a dust protection cover\* over the microscope and store it in a closed container at a dry and mould-free place. Store the microscope and the accessories in the relevant containers when they are not used for a longer time. It is recommended to store also objectives and eyepieces in closed containers with desiccant.

(\*depending on the model a dust protection cover and box can be part of the content)

### NOTICE

**Remember:**

A well maintained microscope will keep its optical quality for years and thus maintain its value.

## 20 Disposal

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

**Do not dispose of electronic devices in the household garbage!**

As per Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

In accordance with the regulations concerning batteries and rechargeable batteries, disposing of them in the normal household waste is explicitly forbidden. Please make sure to dispose of your used batteries as required by law — at a local collection point or in the retail market. Disposal in domestic waste violates the Battery Directive.

Batteries that contain toxins are marked with a sign and a chemical symbol. "Cd" = cadmium, "Hg" = mercury, "Pb" = lead.

## ⚠ CAUTION

### Danger of chemical burns!

Leaking battery acid can lead to chemical burns. Avoid contact of battery acid with skin, eyes and mucous membranes. In the event of contact, rinse the affected region immediately with a plenty of water and seek medical attention.

- Drain the battery completely before disposal.
- Remove the screws in the housing and open the device housing.
- Remove the connections to the battery and remove the battery.
- Cover the open contacts with tape and package the battery securely for transport.

## 21 Technical data

	<b>Biorit TP</b>
<b>Article No.</b>	5101100
<b>Microscope head</b>	monocular
<b>Cross table</b>	Yes, with Nonius scale
<b>Eyepiece(s)</b>	1 x DIN WF 10x
<b>Objective(s)</b>	4x / 10x / 40x
<b>Magnification</b>	40x / 100x / 400x
<b>Power supply</b>	DC 5V / 100 mA
<b>Illumination</b>	LED*

\*The LED illumination used in this device has a very long lifespan and do not need to be changed. LEDs can not be changed due to technical reasons!

## 22 Warranty

The regular guarantee period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required.

You can consult the full guarantee terms as well as information on extending the guarantee period and details of our services at [www.bresser.de/warranty\\_terms](http://www.bresser.de/warranty_terms).

## 23 EC Declaration of Conformity

Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards.

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## GB IE

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail.

**e-mail:** [service@bresseruk.com](mailto:service@bresseruk.com)  
**Telephone\*:** +44 1342 837 098

BRESSER UK Ltd  
Unit 1 starborough Farm,  
Starborough Road, Nr Marsh Green,  
Edenbridge, Kent TN8 5RB  
Great Britain

\*Number charged at local rates in the UK (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

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Folux B.V.